

Technical Bulletin



Anderson Instrument Co.
156 Auriesville Road
Fultonville, NY 12072
1-518-922-5315
Fax 1-800-726-6733
www.andinst.com

Form AIC5024 © 1/04
Effective: July 17, 2006
Supercedes: January 18, 2005

AUTOCLAVEABLE "HA" PRESSURE TRANSMITTERS

- Application:** Sanitary pressure transmitter for autoclaveable equipment/skids.
- Issue:** Sanitary pressure transmitters for autoclave applications are not available.
- Solution:** Following extensive autoclave testing of all models of Anderson "Pharmaceutical Series" pressure transmitters, a modified version of our "HA" pressure transmitter is now available for applications where the pressure sensor must be autoclaved.
- How to Specify:** This transmitter can be specified and priced using the attached matrix. Operating specifications are largely unchanged, with the following provisions:

Process Connections:	1.5" Tri-Clamp® or Anderson "CPM" only
Maximum Autoclave Temperature:	124°C (255°F)
Maximum Autoclave Cycle Time:	1 hour at Max Temp.
Number of Autoclave Cycles:	Warranted for 250 cycles minimum
Cabling:	Quick Disconnect Receptacle (QDR) w/ Field Wireable Connector (FWC) included *

*FWC and attached cabling must be disconnected during autoclave cycles; QDR must be left in place during autoclave cycles.

Please contact our Customer Service Department at 1-800-833-0081 if you have any questions regarding this product. For additional copies of this bulletin, visit our website at www.andinst.com.

